

# TOKU-E Nitrocefin



## Beta-lactamase Detection Protocol

### Principle and Summary

The chromogenic properties of Nitrocefin (TOKU-E item # N005) degradation can be exploited in several simple and rapid assays for detection of  $\beta$ -lactamases. This protocol will outline preparation of Nitrocefin solution and its use in  $\beta$ -lactamase detection methods.

### Solution Preparation

Working solution concentrations of Nitrocefin typically lie within a range of 0.5 mg/ml – 1.0 mg/ml and appear as a yellow color after preparation.

\* Before preparing a working solution, we recommend preparing a concentrated stock solution – approximately 10 mg/ml - in DMSO.

The proceeding example will describe how to prepare a stock solution (10 mg/ml) and a working concentration (1.0 mg/ml) using 10.0 mg of Nitrocefin.

#### Stock solution – 10.0 mg/ml in DMSO

- a. Add 10.0 mg Nitrocefin powder to 1.0 ml DMSO in appropriate container. (If the container only has 10.0 mg Nitrocefin, it may be more efficient to add the DMSO directly to the container, vortex the contents, and add contents to an appropriate container.)

#### Working solution – 1.0 mg/ml in PBS

- a. Add 9.0 mL PBS buffer (100 mM, neutral pH) to the prepared stock solution to bring the final volume to 10.0 ml.

\*If the solution at this concentration produces a red color before exposure to  $\beta$ -lactamase, dilute the working solution 10X or more until a yellow color is achieved.

-Store any remaining stock solution or working solution at -20°C and protect from light. Working solution should be used within 14 days after preparation.

### Rapid Detection Assays

The following assays should produce a positive result within moments after contact with  $\beta$ -lactamase producing organisms. However, it is recommended to allow 20-30 min. before concluding with a negative result.

#### a. Slide Surface Assay

1. Add one drop of 1.0 mg/ml Nitrocefin to the surface of a clean glass slide.
2. Select a colony from an agar surface using a sterile loop and mix with the drop.
3. Appearance of red color within 20-30 min. indicates  $\beta$ -lactamase activity.  
(If necessary, add additional Nitrocefin solution to prevent drying).

#### b. Direct Contact Assay

1. Place one drop of 1.0 mg/ml Nitrocefin directly on the surface of an isolated colony.
2. Appearance of red color within 20-30 min. indicates  $\beta$ -lactamase activity.

c. Broth Suspension Assay


1. Add 3-5 drops of 1.0 mg/ml Nitrocefim to 1 ml of broth suspension.
2. Appearance of red color within 20-30 min. indicates  $\beta$ -lactamase activity

d. Lysed Cell Assay

1. Lyse 1ml of cell suspension by sonication.
2. Add 3-5 drops of 1.0 mg/ml Nitrocefim to lysed cell suspension.
3. Appearance of red color within 20-30 min. indicates  $\beta$ -lactamase activity.

e. Filter Paper Assay

1. Place a small piece of filter paper (~3 x 3 cm) in a clean petri dish or another clean isolated surface and saturate (3-5 ml) with 1.0 mg/ml Nitrocefim
2. Select an isolated colony and smear over the surface of the impregnated filter paper.
3. Appearance of red color within 20-30 min. indicates  $\beta$ -lactamase activity.

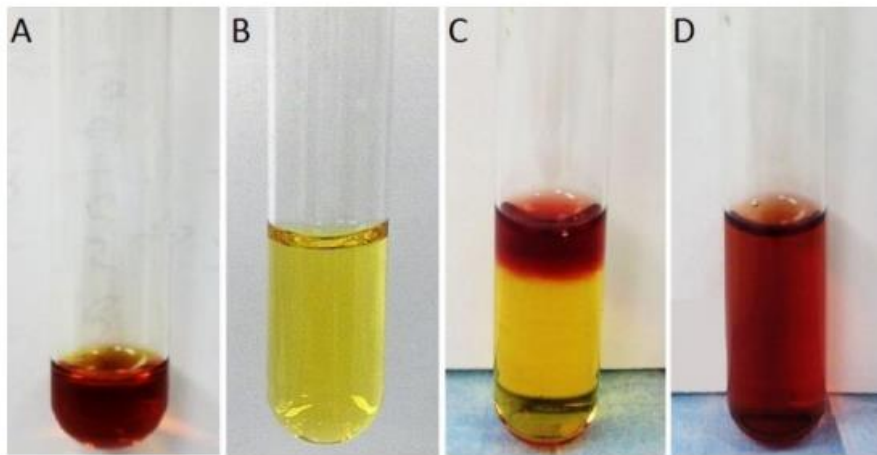

**Fig. 1** Solution preparation and color change before and after  $\beta$ -lactamase exposure

(A) Concentrated nitrocefim (10.0 mg/ml) in DMSO before dilution with PBS buffer.

(B) Nitrocefim diluted with PBS buffer to working concentration (1.0 mg/ml). The yellow color is indicative of intact, undegraded Nitrocefim.

(C) 25 units of  $\beta$ -lactamase dropped on top of Nitrocefim (1.0 mg/mL in PBS). The red color is the result of  $\beta$ -lactamase mediated cleavage of the Nitrocefim.

(D) Vortexed mixture of contents shown in image (C).

### Spectrophotometric analysis of Nitrocefim degradation

In aqueous solution, at approximately neutral pH, Nitrocefim produces two absorption peaks: 217 nm (corresponds to the 7-acyl group) and 386 nm. After exposure to  $\beta$ -lactamase, the peak at 386 nm diminishes, and a new peak appears at 482 nm which corresponds to degraded Nitrocefim. This shift in absorption can be measured over time which enables measurement of the activity of a given  $\beta$ -lactamase enzyme based on the rate at which it degrades Nitrocefim.

The following equations were proposed by O'Callaghan et al.

$$1. \text{ Enzyme activity (pH 7.0, 37}^\circ\text{C, 386nm)} = (x/1.03) \cdot y \cdot V$$

$$2. \text{ Enzyme activity (pH 7.0, 37}^\circ\text{C, 482nm)} = (x/1.59) \cdot y \cdot V$$

Activity is defined as  $\mu\text{mol}$  substrate destroyed per minute per ml of enzyme

$X = \Delta\text{OD}/\text{minute}$

$Y = \text{number of } \mu\text{mol Nitrocefin}$

$V = \text{rate at which Nitrocefin is degraded by a given enzyme (Michaelis constants (Km) and substrate specificity are determined by Line-weaver-Burk plots.}$

1.03 in Equation 1 is equal to the  $\Delta\text{OD}$  (1.75 to 0.72) at 386 nm

1.59 in Equation 2 is equal to the  $\Delta\text{OD}$  (0.15 to 1.74) at 482 nm

## References

O'Callaghan C, Morris A, Kirby SM and Shingler AH (1972) Novel method for detection of  $\beta$  - lactamases by using a chromogenic cephalosporin substrate. *Antimicrob. Agents Chemother.* 1(4):283-288 PMID: [4208895](#)

Parr TR, Pai CH and Bryan LE (1984) Simple screening method for beta-lactamase-positive and-negative ampicillin-resistant *Haemophilus influenzae* isolates. *J. Clin. Microbiol.* 20(1):131-132 PMID [6378964](#)